

NEW

EVO Series, with CMOS detector: StarLine AvaSpec-ULS2048CL-EVO Spectrometer

EVO series: AvaSpec-ULS2048CL



Another new member in our new EVO series: the AvaSpec-ULS2048CL-EVO. Using CMOS technology instead of the conventional CCD technology, this spectrometer offers you the latest technology; ready for the next decade.

The dominant position of CCD detectors in the spectrometer field is fading and new technologies like CMOS have evolved and become a suitable alternative. The AvaSpec-ULS2048CL-EVO offers you this latest technology ensuring a spectrometer platform for the coming years.

In combination with our latest AS-7010 electronics it offers you a versatile device including USB3.0 Communication with 10x higher speed compared to USB2, and a second communication port which offers Gigabit Ethernet for integration in your company network and possibility for long distance communication at an affordable price.

Besides the high speed communica-

tion options, the EVO also offers a fast microprocessor and 50x more memory which can help you to store more spectra onboard and realise more functionality.

Options include a detector collection lens to enhance sensitivity in the 200-1100 nm range and order-sorting filter to reduce 2nd order effects. Furthermore, the AvaSpec-2048CL is available with a wide range of slit sizes, gratings and fiber-optic entrance connectors.

It comes complete with AvaSoft-Basic software, USB cable and an extensive manual.

The AvaSpec-ULS2048CL-EVO is also available as OEM unit, Bench only or Rackmount version.

Technical Data

Optical Bench	ULS Symmetrical Czerny-Turner, 75 mm focal length
Wavelength range	200-1100 nm
Resolution	0.06 -20 nm, depending on configuration (see table)
Stray-light	0.19-1.0%, depending on the grating
Sensitivity	375,000 counts/ μ W per ms integration time
Detector	CMOS linear Image Sensor
Signal/Noise	300:1
AD converter	16-bit, 6 MHz
Integration time	30 μ s - 59s
Interface	USB 3.0 high-speed, 5 Gbps Gigabit Ethernet 1 Gbps
Sample speed with on-board averaging	0.38 ms /scan
Data transfer speed	0.38 ms/scan (USB3), 1.0 ms (ETH)
Digital IO	HD-26 connector, 2 Analog in, 2 Analog out, 13 Digital bidirectional, trigger, sync., strobe, laser
Power supply	Default USB3 power, 500 mA Or 12VDC, 300 mA
Dimensions, weight	177 x 127 x 44,5 mm (1 channel), 1135 grams

Ordering Information

AvaSpec-ULS2048CL-EVO	<ul style="list-style-type: none"> Fiber-optic Spectrometer, 75 mm AvaBench, 2048 pixel CMOS detector 14 x 200 μm, USB powered, high-speed USB 3.0 and ETH interface, incl. AvaSoft-Basic, USB interface cable. Specify grating, wavelength range and options.
PS-12V/1.0A	<ul style="list-style-type: none"> External power supply, needed for operation in ETH mode

Grating selection table for AvaSpec-ULS2048CL-EVO

Use	Useable range (nm)	Spectral range (nm)	Lines/mm	Blaze (nm)	Order code
UV/VIS/NIR	200-1100**	891**	300	300	UA
UV/VIS/NIR	200-1100**	891**	300	300/1000	UNA-DB
UV/VIS	200-850	515	600	300	UB
UV	200-750	247-218*	1200	250	UC
UV	200-650	163-143*	1800	UV	UD
UV	200-580	113-69*	2400	UV	UE
UV	200-400	69-45*	3600	UV	UF
UV/VIS	250-850	515	600	400	BB
VIS/NIR	300-1100**	792**	300	500	VA
VIS	360-1000	495	600	500	VB
VIS	300-800	247-218*	1200	500	VC
VIS	350-750	142-89*	1800	500	VD
VIS	350-640	74-49*	2400	VIS	VE
NIR	500-1050	495	600	750	NB
NIR	500-1050	218-148*	1200	750	NC
NIR	600-1160	346-297	830	800	SI
NIR	600-1100**	495**	300	1000	IA
NIR	600-1100	495	600	1000	IB

* depends on the starting wavelength of the grating; the higher the wavelength, the bigger the dispersion and the smaller the range to select.

** please note that not all 2048 pixels will be used for the useable range

Resolution table (FWHM in nm) for AvaSpec-ULS2048CL-EVO

Grating (lines/mm)	Slit size (μm)					
	10	25	50	100	200	500
300	1.0	1.4	2.5	4.8	9.2	21.3
600	0.40-0.53*	0.7	1.2	2.4	4.6	10.8
830	0.32	0.48	0.93	1.7	3.4	8.5
1200	0.20-0.28*	0.27-0.38*	0.52-0.66*	1.1	2.3	5.4
1800	0.10-0.18*	0.20-0.29*	0.34-0.42*	0.8	1.6	3.6
2400	0.09-0.13*	0.13-0.17*	0.26-0.34*	0.44-0.64*	1.1	2.7
3600	0.06-0.08*	0.10	0.19	0.4	0.8	1.8

* depends on the starting wavelength of the grating; the higher the wavelength, the bigger the dispersion and the better the resolution

Options

-RS	• Replaceable slit
DCL-UV/VIS-200	• Quartz Detector Collection Lens (200-1100 nm)
SLIT-XX	• Slit size, please specify XX = 5, 10, 25, 50, 100, 200 or 500 μm
SLIT-XX-RS	• Replaceable slit with SMA connector, specify slit size XX=25, 50, 100, 200 or 500 μm. Only in combination with AvaSpec-ULS2048CL-EVO-RS
SLIT-XX-RS-FCPC	• as SLIT-XX-RS, but with FC/PC connector
OSF-YYY	• Order-sorting filter for reduction of 2nd order effects please specify YYY= 305, 395, 475, 515, 550 or 600 nm
OSC	• Order-sorting coating with 600 nm long-pass filter for BB (>305 nm) and VB gratings in AvaSpec-ULS2048CL-EVO, recommended with OSF-305
OSC-UA	• Order-sorting coating with 350 and 600 nm long-pass filter for UA, VA gratings in AvaSpec-ULS2048CL-EVO
OSC-UB	• Order-sorting coating with 350 and 600 nm long-pass filter for UB or BB (<350 nm) gratings in AvaSpec-ULS2048CL-EVO
-FCPC	• FC/PC fiber-optic connector