

Levante IR

Tunable OPO for IR Pump Laser

- Levante IR is a synchronously pumped Optical Parametric Oscillator with an innovative and user-friendly design. Mode-locked lasers with an output wavelength of 1 μm are suitable as pump lasers. The output wavelength ranges from approx. 1300 nm to 5000 nm.
- The tunable IR OPO is available as picosecond and femtosecond version. Its synchronous pumping scheme provides a jitter-free and stable repetition rate that is identical to the repetition rate of the pump laser.



- Automated wavelength tuning
- Synchronous pumping scheme
- Standard Software Interface (using TCP/IP)
- Integrated spectrometer for OPO Signal wavelength range

Automated Tuning for IR Wavelength Range

- The output wavelength range for the Signal beam is 1315 ... 2000 nm and the wavelength for the Idler output ranges from 2150 ... 5000 nm. This is ideal for applications that require tunable radiation in the IR range, e.g. vibrational spectroscopy. The wavelength tuning is fully automated.
- An extended wavelength range from 660 nm to 15 μm can be generated when Levante IR is operated in combination with APE's HarmoniXX series. Wavelength converters SHG (Second Harmonic Generation), THG (Third Harmonic Generation) and DFG (Difference Frequency Generation) are available.

Control Software and TCP/IP Standard Software Interface

- All models come with an easy to use control software, allowing for real-time data display. Furthermore, the TCP/IP-based standard software interface by APE makes it straight forward to set up remote control. This allows you, for example, to design your own automated measurement routines. Simply use our protocol templates for rapid configuration with familiar programming languages, including C++, C#, LabVIEW, Python, Matlab, and Ruby.

Software-Interface

- The Levante IR comes with an easy-to-use control interface with real-time data display.



Levante IR Specifications

ps Version

Required Pump laser (not included)	IR ps-Laser 1030 ... 1064 nm 1 ... 10 ps 75 ... > 80 MHz repetition rate (fixed at the time of order)
Wavelength Output (Signal)	1320 ... 2000 nm
Wavelength Output (Idler)	2150 ... 4800 nm (typ. ... 5000 nm)
Power Output at 1500 nm (Signal)	> 1.5 W when pumped with 7.5 W, 2 ps > 2 W when pumped with 10 W, 6 ps
Power Output at 2500 nm (Idler)	> 0.6 W when pumped with 7.5 W, 2 ps > 0.8 W when pumped with 10 W, 6 ps
Pulse Width	typ. 2 ps at 2 ps pump
Time Bandwidth Product	typ. 0.6
Repetition Rate	Equal to pump laser
Wavelength Tuning	Fully automated
Wavelength Tuning Feedback	Built-in spectrometer
Software	Included
Remote Control	Available via TCP/IP

fs Version


Required Pump laser (not included)	IR fs-Laser 1030 ... 1064 nm 70 ... 300 fs 40 ... > 80 MHz repetition rate (fixed at the time of order)
Wavelength Output (Signal)	1320 ... 2000 nm
Wavelength Output (Idler)	2150 ... 4800 nm (typ. ... 5000 nm)
Power Output at 1500 nm (Signal)	> 1.2 W when pumped with 6 W at 1035 nm
Power Output at 2500 nm (Idler)	> 0.6 W when pumped with 6 W at 1035 nm
Pulse Width	typ. 200 fs at 200 fs pump
Time Bandwidth Product	typ. 0.6
Repetition Rate	Equal to pump laser
Wavelength Tuning	Fully automated
Wavelength Tuning Feedback	In-build spectrometer
Software	Included
Remote Control	Available via TCP/IP

... Specifications


Typical Data



Typical Signal power data over Signal wavelength
(pumped with 6 W, ps-Version)

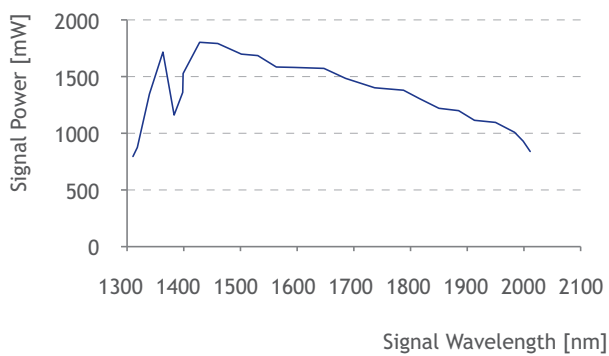

Typical Idler power data over Idler wavelength
(pumped with 6 W, ps-Version)

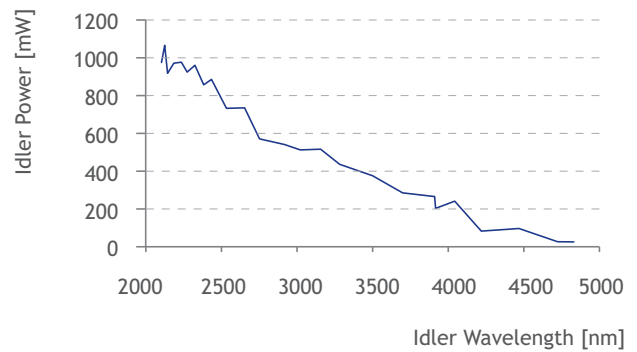

Options

Harmonic Generation SHG or THG for Signal

Harmonic Generation SHG or THG for Idler

Difference Frequency Generation DFG up to 15 μm

Adaption to other pump parameters

Dimensions

1169 mm x 204 mm x 402 mm


Drawings

See appendix for details

Appendix Technical Drawings

All dimensions in mm

Levante IR

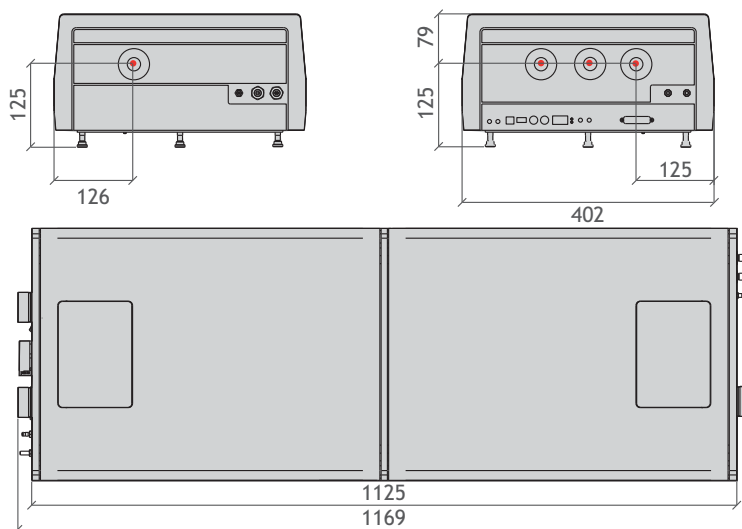


www.phototechnica.co.jp
フォトテクニカ株式会社
 〒336-0017 埼玉県さいたま市南区南浦和 1-2-17
 TEL:048-871-0067 FAX:048-871-0068
 e-mail:voc@phototechnica.co.jp

Contact

APE Angewandte Physik & Elektronik GmbH

Plauener Str. 163-165|Haus N|13053 Berlin|Germany
 T: +49 30 986 011-30
 F: +49 30 986 011-333
 E: sales@ape-berlin.de
 www.ape-berlin.de

APE follows a policy of continued product improvement. Therefore, specifications are subject to change without notice.

© APE GmbH | January 2019 | Rev. 3.1.1