# **ORPHEUS** | PS

## Narrow-Bandwidth Optical Parametric Amplifier

#### **FEATURES**


- Picosecond pulses from femtosecond pump
- 210 4800 nm tuning range
- 800 fs 3 ps pulse duration
- < 20 cm<sup>-1</sup> spectral bandwidth
- Up to 100 kHz repetition rate
- High output stability



ORPHEUS-PS is a narrow-bandwidth optical parametric amplifier (OPA), designed for PHAROS and CARBIDE lasers. The device is pumped by picosecond pulses produced in a second harmonic bandwidth compressor SHBC and is seeded by a white-light continuum generated by femtosecond pulses. This enables very high pulse-to-pulse energy stability compared to other methods for generating tunable picosecond pulses. The white-light generation is integrated into the same housing as the amplification, enabling high long-term stability and

ease of use. The system features high conversion efficiency, bandwidth- and diffraction-limited output, and full computer control.

Part of the pump laser radiation can be split to simultaneously pump a femtosecond OPA, giving access to the complete set of beams necessary for versatile spectroscopy applications such as femtosecond stimulated Raman spectroscopy (FSRS) and sum-frequency generation (SFG) spectroscopy.

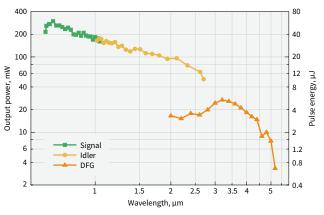

**ORPHEUS-PS** tuning curves. Pump: 5 W, 1000 μJ, 5 kHz from PHAROS-SP.




#### **SPECIFICATIONS**


Model	ORPHEUS-PS
MAIN OUTPUT	
Tuning range 1)	640 – 1000 nm (Signal) 1060 – 2600 nm (Idler)
Conversion efficiency at peak	> 8% (Signal and Idler combined)
Pulse duration	800 fs – 3 ps
Spectral bandwidth	< 20 cm <sup>-1</sup> @ 700 – 2000 nm
Pulse-to-pulse energy stability <sup>2)</sup>	< 2% @ 800 nm
AUXILIARY OUTPUT 1 (515 nm)	
Center wavelength 3)	515 nm ± 5 nm
Generation efficiency 4)	> 15%
AUXILIARY OUTPUT 2 (1030 nm)  Center wavelength 5)	1030 ± 10 nm
Pulse duration	< 300 fs
Pulse energy	> 5 μJ
WAVELENGTH EXTENSION	
SH package at peak (320 – 500 nm (SHS), 530 – 640 nm (SHI))	> 3%
FH package at peak (210 – 250 nm (FHS), 265 – 320 nm (FHI))	> 0.3% <sup>6)</sup>
2400 – 4800 nm (DFG)	> 0.25% @ 3200 nm <sup>6)</sup>
PUMP LASER REQUIREMENTS	
Pump source	PHAROS or CARBIDE with uncompressed output option
Wavelength	1030 nm
Repetition rate	Single-shot – 100 kHz
Maximum pump power	20 W
Pump pulse energy	100 μJ – 3.2 mJ

- <sup>2)</sup> Expressed as NRMSD (normalized root mean squared deviation).
- <sup>3)</sup> Direct SHBC output, not simultaneous to OPA; see details in SHBC
- 5) Compressed pump output.
- <sup>6)</sup> For > 200 μJ pump pulse energy.



### **DRAWINGS**

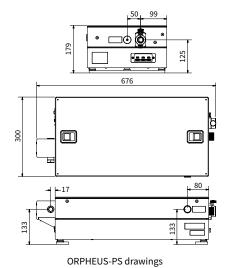

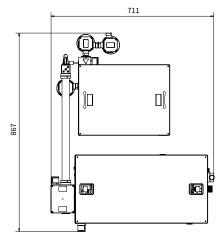

ORPHEUS-PS with SHBC drawing



〒336-0017 埼玉県さいたま市南区南浦和 1-2-17 TEL:048-871-0067 FAX:048-871-0068 e-mail:voc@phototechnica.co.jp

